

Room Type/Surface	Minimum Frequency* More frequent cleaning may be required depending on conditions	Method	Application	Product (must be sterile)
ISO 5 PRIMARY ENGINEERING CONTROLS				
<ul style="list-style-type: none"> (LAFW) Laminar Air Flow Workstations (BSC) Biological Safety Cabinet (CAI) Compounding Aseptic Isolator (CACI) Compounding Aseptic Containment Isolator 	<ul style="list-style-type: none"> Beginning of each shift 	Wipe	Cleaning/Disinfecting	Step 1. Decon-Quat 200C Step 2. Decon-Ahol 70% IPA <i>or</i> Alcoh-Wipe 70% IPA
	<ul style="list-style-type: none"> Before each batch Every 30 minutes during compounding When spills occur When surface contamination is known/suspected 	Wipe	Disinfecting	Decon-Ahol 70% IPA <i>or</i> Alcoh-Wipe 70% IPA
	<ul style="list-style-type: none"> Monthly or quarterly based in response to EM micro monitoring results 	Wipe	Cleaning/Sporicidal	Step 1. Decon-Clean spray <i>or</i> wipe Step 2. Decon-Spore 200 Plus <i>or</i> Hypo-Chlor 5.25% <i>or</i> Hypo-Chlor 0.52% <i>or</i> Steri-Perox 6% spray <i>or</i> wipe
	<ul style="list-style-type: none"> During use as required 	Wipe	Cleaning	Decon-Ahol 70% IPA
ISO 7 & 8 BUFFER AND ANTEROOM				
<ul style="list-style-type: none"> Counters Work surfaces Door plates Handles Gowning benches 	<ul style="list-style-type: none"> Daily 	Wipe	Cleaning/Disinfecting	Decon-Quat 200C
<ul style="list-style-type: none"> Floors 	<ul style="list-style-type: none"> Daily (when no aseptic operations are processing) 	Mop	Cleaning/Disinfecting	Decon-Quat 200C
	<ul style="list-style-type: none"> Monthly or quarterly based in response to EM micro monitoring results 	Mop	Cleaning/Sporicidal	Step 1. Decon-Clean spray <i>or</i> wipe Step 2. Decon-Spore 200 Plus <i>or</i> Hypo-Chlor 5.25% <i>or</i> Hypo-Chlor 0.52% <i>or</i> Steri-Perox 6% spray <i>or</i> wipe
<ul style="list-style-type: none"> Walls and Ceilings 	<ul style="list-style-type: none"> Monthly 	Mop	Cleaning/Disinfecting	Decon-Quat 200C
	<ul style="list-style-type: none"> Monthly or quarterly based in response to EM micro monitoring results 	Mop	Cleaning/Sporicidal	Step 1. Decon-Clean spray <i>or</i> wipe Step 2. Decon-Spore 200 Plus <i>or</i> Hypo-Chlor 5.25% <i>or</i> Hypo-Chlor 0.52% <i>or</i> Steri-Perox 6% spray <i>or</i> wipe
<ul style="list-style-type: none"> Storage Shelving 	<ul style="list-style-type: none"> Monthly 	Wipe	Cleaning/Disinfecting	Decon-Quat 200C
<ul style="list-style-type: none"> Incoming Supplies 	<ul style="list-style-type: none"> As needed 	Wipe	Disinfecting	Step 1. Decon-Quat 200C Step 2. Decon-Ahol 70% IPA <i>or</i> Alcoh-Wipe 70% IPA
CHEMO COMPOUNDING				
<ul style="list-style-type: none"> Deactivate Chemo Drugs 	<ul style="list-style-type: none"> As needed 	Wipe	Deactivate/Disinfecting/Rinsing	Step 1. Hypo-Chlor 5.25% Step 2. Decon-Clean spray <i>or</i> wipe Step 3. Steri-Water <i>or</i> Decon-Ahol 70% IPA <i>or</i> Alcoh-Wipe 70% IPA

Always refer to site specific SOP's for cleaning and disinfection instructions.

Product	Dilution	Application
Decon-Quat 200C #19178, 19179, 19180	1:128 RTU	Cleaning/Disinfecting
Decon-Ahol 70% IPA #19175, 19176	RTU	Disinfecting
Alcoh-Wipe 70% IPA #19174	RTU	Disinfecting
Decon-Clean #19177,19333	1:128 RTU	Cleaning
Decon-Spore 200 Plus #19181	6.4:128 RTU	Sporicidal/Disinfecting
Hypo-Chlor 5.25% spray <i>or</i> wipe #19183, 18733	0.52% RTU	Cleaning/Disinfecting
Hypo-Chlor 0.52% spray <i>or</i> wipe #19183, 19184	0.52% RTU	Cleaning/Disinfecting
Steri-Perox 6% spray <i>or</i> wipe #19188, 19189	RTU	Decontamination
Wipes #10091, 18956, 18957	N/A	Wiping

NOTE: All products used in the critical classified area must be sterile. Product information by Veltek Associates, Inc.