



VELTEK ASSOCIATES, INC.

# TECHNICAL DATA FILES



## **HYPO-CHLOR<sup>®</sup> Wipe 0.25% & 0.52%, and 5.25%**

Sodium Hypochlorite and Water for Injection Solution  
Pre-Saturated 100% Polyester Wipe

## Product Overview

**HYPO-CHLOR® Wipes** are saturated with a ready-to-use 0.25%, 0.52%, or 5.25% concentration of sodium hypochlorite solution formulated with USP Water for Injection. These wipes can be used in both aseptic and non-aseptic wipe downs of filling and packaging machinery, stainless steel surfaces, lexan, polycarbonate, glass, and any critical surface that requires cleaning. Ideal use is in broad or intricate locations where a sodium hypochlorite spray could create an excess of residue. Pre-saturated wipers, including **HYPO-CHLOR Wipes**, can prevent the overuse of chemicals and reduce costs. **HYPO-CHLOR Wipes** are available in a standard 12”x12” (30.5 cm x 30.5 cm) size and either sterile or non-sterile.

**HYPO-CHLOR Wipes** are manufactured via aseptic fill at 0.2 microns into gamma irradiated sterile components in ISO 5 (Grade A/B, Former Class 100). Individual lots are lot sterility tested according to current USP compendium, are completely traceable, and completely validated for sterility and shelf life. **HYPO-CHLOR Wipes** are delivered each time with lot specific Certificate of Analysis, Certificate of Sterility, and Certificate of Conformance tested to current USP compendium.

The material used in our **HYPO-CHLOR Wipes** is a 100% continuous knit polyester filament that has been specifically designed to be exceptionally clean and absorbent. Particulates, residues, and foreign matter is effectively entrapped in the webbing of the material. The knit construction and pattern has been developed to ensure the greatest product integrity and cleaning performance while offering high resistance to abrasion and picking.

FocusEdge cutting technology is the standard that VAI® uses to cut their wipes. This cutting method does not heat the yarn bundles as high as lasers, therefore, eliminating the small plastic balls that form on the end of every yarn strand. Due to the wipe’s sealed edges, shedding is minimal during wipe downs or general use, therefore, reducing particulates in controlled environments. After cutting, the wipes are then laundered in ISO Class 4 conditions to ensure cleanliness.

VAI’s **HYPO-CHLOR Wipes** are packaged in an opaque, zip-lock style, ported bag that allows for continued use after opening. Each sterile wiper bag is double bag packaged and available in our ABCD Cleanroom Introduction System® for easy introduction into the core or other critical environments. In addition, **HYPO-CHLOR Wipe 0.52%** is available individually packaged in an opaque, easy-tear style bag.

## Product Uses

- Aseptic or non-aseptic environments
- Wipe downs on filling and packing machinery
- Spill control
- On lexan, glass, polycarbonate, stainless steel or any other critical surface that requires cleaning
- Where a sodium hypochlorite spray would create excess residue
- To prevent overuse of sodium hypochlorite when compared to using trigger sprayers or gallon pour



## Quality and Manufacturing

- Wipe material laundered in ISO 4 (Grade A/B, Former Class 10)
- Sodium hypochlorite solution is filtered at 0.2 microns
- Formulated with USP Water for Injection (0.25 EU/mL)
- Filled in ISO 5 (Grade A/B, Former Class 100)
- Cut using FocusEdge cutting technology
- Aseptically filled into sterile components via gamma irradiation at  $10^{-6}$  SAL
- Lot sterility tested according to current USP Compendium
- Completely lot traceable
- Completely validated for sterility and shelf life



Ported bag for aseptic filling into sterile components

## Features and Benefits

- ISO 5 (Grade A/B, Former Class 100) Cleanliness
- Each sterile bag is double bagged in easy tear bags
- Quadruple bagged via the ABCD Cleanroom Introduction System®
- Packaged in an opaque, zip-lock style, easy tear bag
- Re-sealable for continued use after opening
- Individually labeled with lot number and expiration
- Three concentrations: 0.25%, 0.52%, or 5.25% to fit your specific needs
- For use on multiple surfaces and ready-to-use
- Wipes folded for easy removal
- 100% synthetic filament polyester constructed in a continuous knit that minimizes fiber and particulate release
- Material is exceptionally clean and absorbent
- Resistance to abrasion and picking
- FocusEdge allows for minimal shedding via sealed edges
- Standard 12”x12” (30.5 cm x 30.5 cm) size
- Delivered with lot specific Certificate of Analysis, Certificate of Sterility, and Certificate of Conformance tested to current USP Compendium



Easy Tear Packaging



Zip-Lock Style Bag

## ABCD Cleanroom Introduction System®

The ABCD Cleanroom Introduction System is a packaging system that allows operators/users to take the package through each level of classified areas by simply removing one bag at a time. Each bag acts as barrier protecting the finished product from becoming a carrier of viable and non-viable contamination. This prevents the need to decontaminate each outer bag prior to entering a cleaner area. In this packaging system, sterilized groups of containers are contained in two outer bags and after each are removed individual containers are each additionally contained in two easy tear bags.

## Material Specifications

| HYPO-CHLOR Wipe – Laundered Polyester Wipe |   |
|--|---|
| Material Details                           | Specification                                 |
| Material:                                  | 100% Polyester                                |
| Material Description:                      | Laundered Polyester Wipe                      |
| Color:                                     | White   |
| Classification Compatibility:              | ISO Class 5 (Former Class 100)                |
| Construction:                              | Continuous Filament Knitted Construction      |
| Dimension:                                 | 12”x12” (30.5 cm x 30.5 cm)                   |
| Thickness:                                 | 0.56 – 0.64 mm                                |
| Sorption Capacity:                         | >/= 270 mL/m <sup>2</sup>                     |
| Particles:                                 | <4.5x10 <sup>6</sup> particles/m <sup>2</sup> |
| Fibers @ 100 um/m <sup>2</sup> :           | 63, 1.0 IEST-RP-CC004.3, SEC 6.1.3/6.2.2      |

\*Treatable surface area not to exceed 1 sq. meter



100% Polyester Wiper Substrate

## Chemical Specifications

| HYPO-CHLOR Wipe – Sodium Hypochlorite and Water for Injection |                   |
|---|-------------------|
| Certificate of Analysis                                       | Specifications    |
| Assay 0.25%:  | 0.25% - 0.30% w/w |
| Assay 0.52%:  | 0.4% – 0.6% w/w   |
| Assay 5.25%   | 4.0% - 6.0% w/w   |
| Litmus Paper Turns Blue:                                      | Pass              |
| Addition of HCL Gives Off CL <sub>2</sub> Gas:                | Pass              |
| Yellow Flame Test:  | Pass              |
| Expiration Period:  | 18 months         |

## Ordering Information

| HYPO-CHLOR Wipe – Sodium Hypochlorite and Water for Injection |  |          |
|---|--|----------|
| Part number   | Description  | Qty/Case |
| VEL9-12X12-S-3020   | HYPO-CHLOR® Wipe 0.25%, 20 wipes/bag, 10 bags/case, Sterile    | 200      |
| VEL9-12X12-3021   | HYPO-CHLOR Wipe 0.25%, 20 wipes/bag, 10 bags/case, Non-Sterile | 200      |
| VEL9-12X12-S-3018   | HYPO-CHLOR Wipe 0.52%, 20 wipes/bag, 10 bags/case, Sterile     | 200      |
| VEL9-12X12-3019   | HYPO-CHLOR Wipe 0.52%, 20 wipes/bag, 10 bags/case, Non-Sterile | 200      |
| VEL9-12X12-S-3026   | HYPO-CHLOR Wipe 0.52%, Individually Packaged, Sterile          | 100      |
| VEL9-12X12-S-3025   | HYPO-CHLOR Wipe 5.25%, 20 wipes/bag, 5 bags/case, Sterile      | 100      |



VEL9-12X12-3021



VEL9-12X12-S-3026



VEL9-12X12-S-3025

# PRODUCT LABELING

## HYPO-CHLOR® Wipe 0.25%, 0.52%, & 5.25% Sodium Hypochlorite and Water for Injection Solution

(Any specific product label is available upon request.)



**HYPO-CHLOR Wipe Family**

Veltek Associates, Inc.

15 Lee Boulevard, Malvern, PA 19355-1234

Tel: 610-644-8335 Fax: 610-644-8335 [www.sterile.com](http://www.sterile.com) Rev: 05Aug2021

## HYPO-CHLOR® Wipe 0.25% and 0.52%

**Sterile Pharmaceutical Cleanroom Wipe 30.5cm x 30.5 cm (12 in x 12 in) Pre-saturated  
Laundered Polyester Wipe**

**Ready to Use Wipe**

### HYPO-CHLOR Wipe 0.25%

**Ingredients 0.25%:**

Sodium Hypochlorite..... (CAS 7681-52-9)  
.....0.25%  
Other Ingredients: Water

Case# 0.25%: VEL9-025-00 • Rev 1/21

Case# 0.52%: VEL9-052-00 • Rev 1/21

Patents: [www.sterile.com/patents](http://www.sterile.com/patents).

Made in Korea to VAI quality standards.

**Manufactured by:**

Veltek Associates, Inc.  
15 Lee Boulevard  
Malvern, PA 19355-1234 USA  
Tel: 1-610-644-8335 • [www.sterile.com](http://www.sterile.com)

### HYPO-CHLOR Wipe 0.52%

**Ingredients 0.52%:**

Sodium Hypochlorite..... (CAS 7681-52-9)  
.....0.52%  
Other Ingredients: Water

Veltek Associates, Inc.  
Postbus 1062. 8200 BB  
Lelystad, Nederland  
Tel: 00-800-00888700 • [www.sterile.com](http://www.sterile.com)

Harmful to aquatic life with long lasting effects.

**PRECAUTIONS FOR USE:**

Avoid release to the environment. Dispose of contents and container to an authorised waste collection point.

IRL PCS#: 98537

See container for lot# / Exp Date

SDS 0.25%: VEL-130 (Americas), VEL-130-EU (Europe)

SDS 0.52%: VEL-131 (Americas), VEL-131-EU (Europe)

**EMERGENCY NUMBER:** 1-866-928-0789 (*toll free USA*), 1-800-579-7421 (*toll free Canada*), 1-215-207-0061 (*Americas*), +44 1235 239670 (*Europe*), +44 1235 239671 (*Middle East/Africa*), +353 (0)1 809 2166 (*Ireland, available to the public 08.00-22.00*), 000 800 100 7479 (*Hindi*).

**See Safety Data Sheet for first aid and additional product information.**

**For Industrial Use Only.**



## HYPO-CHLOR® Wipe 5.25%

Pharmaceutical Cleanroom Wipe, 12in x 12in (30.5cm x 30.5 cm)

Sodium Hypochlorite at 5.25% wt./wt. in USP Water for Injection

### Ingredients 5.25%:

Sodium Hypochlorite (CAS-No 7681-52-9) .....(525 mL).....5.25%



**DANGER**

Case#: VEL9-5.25-00 • Rev 10/19

Patents: [www.sterile.com/patents](http://www.sterile.com/patents).

Made in Korea to VAI quality standards.

### Manufactured by:

Veltek Associates, Inc.  
15 Lee Blvd. Malvern, PA 19355-1234 USA  
Tel: 1-610-644-8335  
[www.sterile.com](http://www.sterile.com)

Veltek Associates, Inc.  
Postbus 1062 8200 BB Lelystad,  
Nederland  
Tel: +800-00888700  
[www.sterile.com](http://www.sterile.com)

**DANGER.** May be corrosive to metals. Causes skin irritation. Causes serious eye damage. Very toxic to aquatic life with long lasting effects.

**PRECAUTIONS FOR USE:** Keep only in original packaging. Wear eye protection, protective gloves. **IF IN EYES:** Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Immediately call a doctor. Collect spillage. Dispose of contents and container to an authorised waste collection point. Contact with acid liberates toxic gas.

SDS: VEL-119 (America), VEL-119-EU (Europe)

**Emergency Number:** 1-866-928-0789 (toll free USA), 1-800-579-7421 (toll free Canada), 1-215-207-0061 (Americas), +44 1235 239670 (Europe), +44 1235 239671 (Middle East/Africa). +353 (0)1 809 2166 (Ireland) (Available to the public 08.00-22.00), 000 800 100 7479 (Hindi).

**See Safety Data Sheet for first aid and additional product information.**

**For Industrial Use Only.**



## Additional Documentation

Upon request, the following additional documentation is available:

- Specific Product Testing Reports
- Safety Data Sheet
  - **HYPO-CHLOR Wipe 0.25%** - SDS# VEL-131
  - **HYPO-CHLOR Wipe 0.52%** - SDS# VEL-130
  - **HYPO-CHLOR Wipe 5.25%** - SDS# VEL-119
- Product Validation
- Sample lot specific documentation packages including Certificates of Sterility, Certificates of Analysis, and Certificates of Conformance



VAI®'s Wiper Family

Veltek Associates, Inc. *Disposable Products Manufacturing Division*'s capabilities include sterile disposable garments, sterile disposable face masks, dry wipers, sterile wipers, custom cleanroom documentation, sterile synthetic writing substrate, and a HEPA filtered cleanroom printer. DPMD has responded and met demands of professionals in the industry since 1985. Our manufacturing capabilities assure that we will be able to meet all requests and fulfill all of our customer's needs. All DPMD products are processed and packaged in cleanroom facilities from beginning to end and subsequently gamma irradiated or aseptically filled in order assure sterility. For more information call 610-644-8335, email [sales@sterile.com](mailto:sales@sterile.com) or visit our website at [www.sterile.com](http://www.sterile.com).

Patents: [www.sterile.com/patents](http://www.sterile.com/patents)