


INSTRUCTIONS FOR VIAL DECAPPER PLIERS (HCL #7713, 7714, 7718-7720)

To remove crimped aluminum seals on glass vials:

- A. Insert vial squarely into open decapper jaws.
- B. Squeeze firmly.
- C. Holding vial upright, tip decapper to remove aluminum seal.

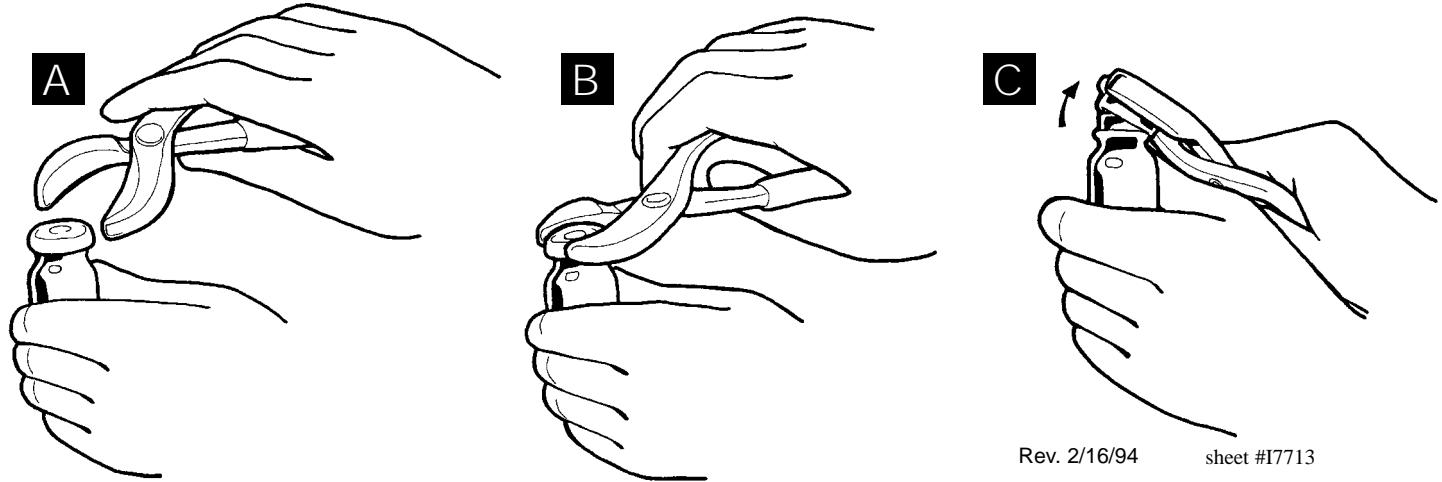

Rev. 2/16/94

sheet #17713

Call Free 1-800-848-1633 • Fax Free 1-800-447-2923
International Call (740) 477-1686 • International Fax (740) 477-2923